

pVectOZ-GFP (Green Fluorescent Protein Expression Vector)

Description

The pVectOZ-GFP vector has been created to produce the highest levels of Green Fluorescent Protein expression in a broad range of mammalian cells and tissues. It contains a proprietary modified human cytomegalovirus (CMV) promoter followed by a specific intron, enhancer and a terminator. The expression vector is engineered in an optimized plasmid backbone to achieve the highest levels of transgene expression in mammalian cells and high copy number production in *Escherichia coli*.

Kit contents

Ref. #PL00120
100µg pVectOZ-GFP (encoding for Green Fluorescent Protein) plasmid in 100µl sterile TE buffer.

Storage

Store at -20°C.

Selection Marker

Kanamycin is the selection gene included for producing the plasmid in *Escherichia coli*.

Applications

pVectOZ-GFP (Green Fluorescent Protein) vector is suitable for all transfection applications (*in vitro* & *in vivo*).

Presentation. The transgene expression level depends mainly on the promoter, enhancer, terminator and plasmid backbone. The pVectOZ-GFP expression cassette was designed to express very high levels of transgene product in many mammalian cells and tissues. This vector has been modified to eliminate sequences affecting transgene expression levels while optimizing those critical for high levels of expression. The final expression cassette accommodates the high levels of transgene expression in mammalian cells as well as high yield of plasmid production in *Escherichia coli*. The

resulting plasmid is the ideal vector to reach high levels of expression *in vitro* and *in vivo*.

Use. For high levels of transgene expression in mammalian cells and tissues. For optimal results, this vector can be used with all OZ Biosciences transfection reagents to transfect a wide variety of mammalian cells and tissues.

GFP detection

For transfections performed with pVectOZ-GFP plasmids, the detection can be done by fluorescent (epifluorescence) or confocal microscopy. The GFP produced has an excitation peak at 470-480 nm and emission peak at 510 nm. GFP expression level can also be monitored by fluorescence-activated cell sorter analysis (FACS).

References

1. Kain, SR. *et al.* (1995). *Biotechniques*. **19**: 650.
2. Cheng, L. *et al.* (1996). *Nature Biotech.* **14**: 606.
3. Welsh, S. *et al.* (1997). *Curr Opin Biotechnol.* **8**: 617.
4. Zelphati, O. *et al.* (1999). *Hum Gene Ther.* **10**: 15.
5. Ward, T. *et al.* (2006). *Meth Biochem Anal.* **47**: 305.

Contact Us

Please, feel free to contact us for all complementary information and remember to visit our website (www.ozbiosciences.com) to stay informed on the latest breakthrough technologies and updated on our complete product list.

Technical questions: tech@ozbiosciences.com
Order: order@ozbiosciences.com
General Information: contact@ozbiosciences.com


OZ Biosciences SAS

163 avenue de Luminy
Case 922, zone entreprise
13288 Marseille cedex 09 -
FRANCE
Ph: +33 (0) 486 948 516
Fax: +33 (0) 486 948 515
contact@ozbiosciences.com
order@ozbiosciences.com

OZ Biosciences INC

4901 Morena Blvd,
Suite 901
San Diego CA 92117 - USA
Ph : + 1-858-246-7840
Fax : + 1-855-631-0626
contactUSA@ozbiosciences.com
orderUSA@ozbiosciences.com

Plasmid Map

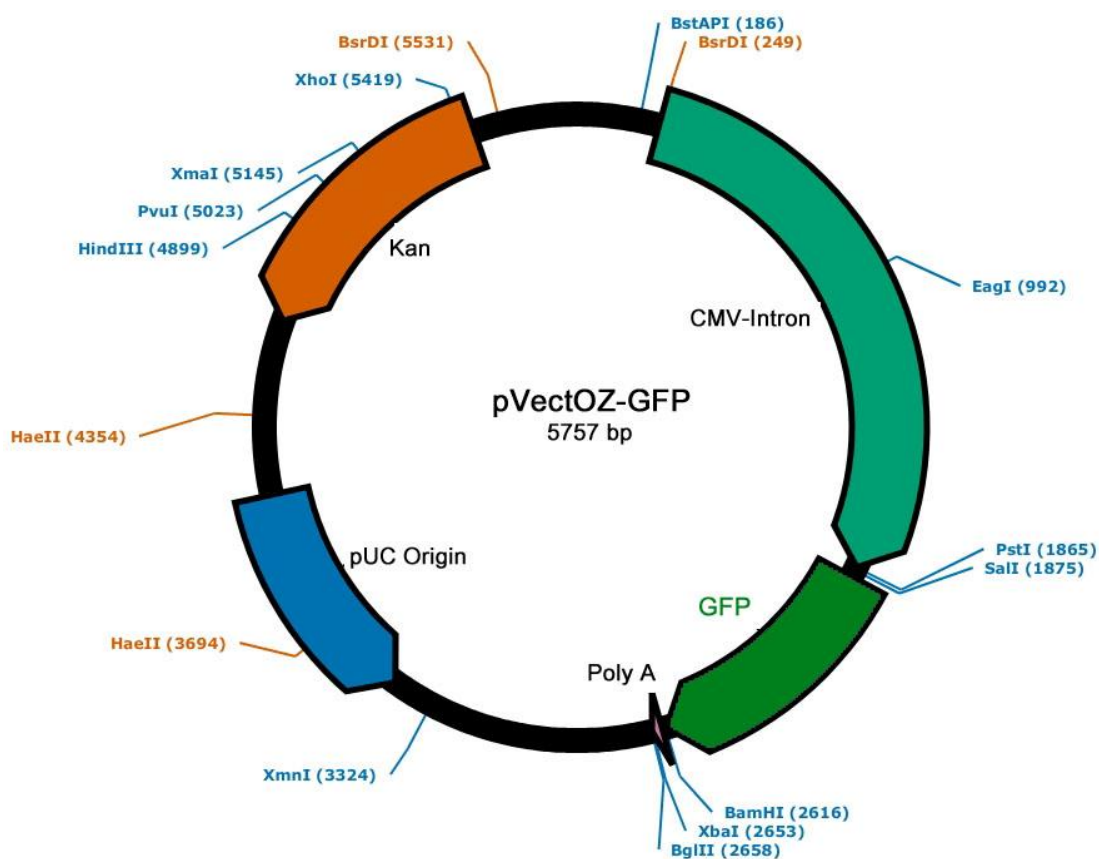

Related Products

Description	Reference	Description	Reference
Magnetofection Technology		Gene & Protein Tools	
Mega Magnetic Plate	MF14000	Bradford – Protein Assay Kit	BA00100
Super Magnetic Plate	MF10000	GeneBlaster selection kit	GB20010
Magnetic Plate 96-magnets	MF10096	β-Galactosidase (ONPG) assay kit	GO10001
PolyMag 1 mL (for all nucleic acids)	PN31000	β-Galactosidase (CPRG) assay kit	GC10002
PolyMag Neo 1 mL (for all nucleic acids)	PG61000	X-Gal Staining Kit	GX10003
CombiMag 1 mL (boost transfection reagent)	CM21000	Plasmids	
SilenceMag 500μL (for siRNA applications)	SM10500	pVectOZ-CAT 25μg	PL00010
NeuroMag 1 mL (for neuron transfection)	NM51000	pVectOZ-GFP 25μg	PL00020
Lipofection (lipid-based reagents)		pVectOZ-LacZ 25μg	PL00030
DreamFect Gold Transfection reagent 1mL	DG81000	pVectOZ-Luc 25μg	PL00040
DreamFect Transfection reagent 1mL	DF41000	pVectOZ-SEAP 25μg	PL00050
Lullaby siRNA Transfection reagent 1mL	LL71000	pVectOZ-CAT 100μg	PL00110
VeroFect Transfection Reagent 1mL	VF61000	pVectOZ-LacZ 100μg	PL00130
FlyFectin Transfection Reagent 1mL	FF51000	pVectOZ-Luc 100μg	PL00140
		pVectOZ-SEAP 100μg	PL00150