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Lysozyme

from hen egg white

Synonyme(s):
Mucopolysaccharide N-acetylmuramoyl-hydrolase, Muramidase, lysozyme

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
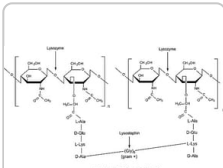
L6876

Lysozyme from chicken egg white

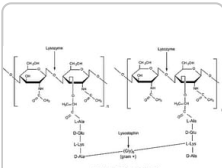lyophilized powder, protein ≥90 %, ≥40,000 units/mg

Sigma-Aldrich

L4919

Lysozyme from chicken egg white

BioUltra, lyophilized powder, ≥98% (SDS-PAGE),

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DESCRIPTION

Description générale

Muramidase, Mucopeptide N-acetylmuramoyl-hydrolase

Application

Lysozyme has been used for:

- Cell wall degradation
- Preparation of protoplasts
- Bacteriolysis
- Pharmacology (anti-inflammatory effect on mucous membrane; healing of tissue; protection against bleeding; normalization of mucus secretion; dissolution of pyrogens; potentiation of antibiotics)
- Food and drinks (flavor enhancer)
- Sample preparation prior to isolation of nucleic acids

Lysozyme has been used in the monitoring of transcription by real-time -PCR and sample preparation.

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Lysozyme hydrolytically cleaves the bonds between N-acetyl-D-glucosamine and N-acetylmuramic acid residues (GlcNAc β 1-4Mur) in mucopolysaccharides and the glycan skeletons of the corresponding peptidoglycans (the latter are glycosaminoglycans linked to or crosslinked by peptide chains). Lysozyme effectively destroys the cell walls of bacteria.

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One Shugar unit is that quantity of enzyme in 1 ml of a suspension of *Micrococcus luteus* of pH 7.0 whose initial absorbance at a wavelength of 450 nm is 0.750 for a pathlength of 10 mm which causes the absorbance to decrease at the rate of 0.001 per minute, all measurements being carried out at a temperature of +25 °C.

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Working solution: Lysozyme hydrochloride is readily soluble in water and buffer solutions but insoluble in organic solvents.

Storage conditions (working solution): -15 to -25 °C

Aqueous solutions of the hydrochloride (2 mg/ml dist. water) can be stored for several days at 2 to 8° C or at -15 to -25° C for several weeks.

Reconstitution

Prepare a stock solution at a concentration of 50 mg/ml in dist. water. Dispense into aliquots and store at -15 to -25 °C. Discard each aliquot after use.

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Store at 2 to 8 °C. (Store dry!)

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For life science research only. Not for use in diagnostic procedures.

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
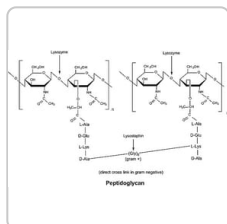
L8402

Lysozyme from human neutrophils

≥95% (SDS-PAGE), lyophilized powder,
≥100,000 units/mg protein (E1%/280)

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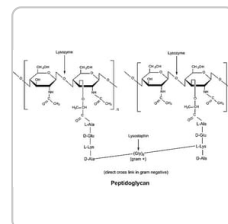Modifications to the peptidoglycan backbone help bacteria to establish infection.

Kimberly M Davis et al.

Infection and immunity, 79(2), 562-570 (2010-11-03)

Bacterial pathogens that colonize mucosal surfaces have acquired resistance to antimicrobials that are abundant at these sites. One of the main antimicrobials present on mucosal surfaces is lysozyme, a muramidase that hydrolyzes the peptidoglycan backbone of bacteria. Cleavage of the

A new iron acquisition system in Bacteroidetes.

Manfredi P, et al.

Infection and Immunity (2014)

Visualization of RNA 3' ends in Escherichia coli Using 3' RACE Combined with Primer Extension

Xun Wang, et al.

Bio-protocol, 8(5) (2018)

A computationally designed chimeric antigen receptor provides a small-molecule safety switch for T-cell therapy.

Greta Giordano-Attianese et al.

Nature biotechnology, 38(4), 426-432 (2020-02-06)

Approaches to increase the activity of chimeric antigen receptor (CAR)-T cells against solid tumors may also increase the risk of toxicity and other side effects. To improve the safety of CAR-T-cell therapy, we computationally designed a chemically disruptable heterodimer (CDH)

Isolation of Heterocysts from *Anabaena* sp. PCC 7120

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Roche
FORMD-RO
Formate Dehydrogenase
from *Candida boidinii*

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.