[Toutes les photos \(2\)](#)

Documents

[↓ FDS](#)[🔍 COO/COA](#)[Plus de documents »](#)

ENDOGLUS-RO Roche

Endoproteinase Glu-C Sequencing Grade

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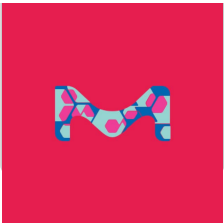
P2922

Endoproteinase Glu-C from *Staphylococcus aureus* V8

Type XVII-B, lyophilized powder, 500-1,000 units/mg solid

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10791156001

Endoproteinase Glu-C (V8 Protease)
from *Staphylococcus aureus* V8

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DESCRIPTION

Description générale

Endoproteinase Glu-C Sequencing Grade is a serine protease isolated from *Staphylococcus aureus* V8 as a highly purified and specific protease. Inhibitors of Endoproteinase Glu-C are DFP, α 2-macroglobulin and TLCK.

Spécificité

Serine protease that very specifically hydrolyzes peptide bonds at the carboxylic side of Glu (in ammonium bicarbonate, pH 7.8, or ammonium acetate buffer, pH 4.0), or Glu and Asp (in phosphate buffer, pH 7.8). The specificity of Endoproteinase Glu-C Sequencing Grade is verified with the oxidized B-chain of insulin (insulin B_{ox}) as substrate. High concentrations of Endoproteinase Glu-C Sequencing Grade (1 part by weight enzyme with 10 parts by weight insulin B_{ox}) are incubated to detect traces of impurities, for example, contaminating proteases.

Heat inactivation: Most proteases sequencing grade are stopped by boiling for 5 minutes.

Application

Use Endoproteinase Glu-C Sequencing Grade for protein structure analysis and for sequence analysis.

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Free of impurities that may interfere with separation of peptides in RP HPLC.

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Working concentration: 1/100 to 1/20 (w/w)

With a minimum amount of 1 to 5 µg insulin B_{ox}, sufficient levels can be detected using the Column Aquapore RUP 300, 4.6 x 100 nm, 7 µm.

Working solution: Solvent is recommended in double-distilled water.

Solvents other than water have not been tested; Roche recommends using 50 mM Hepes, pH 7.8.

Storage conditions (working solution): +2°C to +8°C for 1-2 days.

Stability: At +2 to +8°C within specification range for 12 months; stored dry.

Reconstitution

Lyophilized Endoproteinase Glu-C sequencing grade is reconstituted in double-dist. water. To avoid autolysis, the incubation temperature should not exceed 25 °C.

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For life science research only. Not for use in diagnostic procedures.

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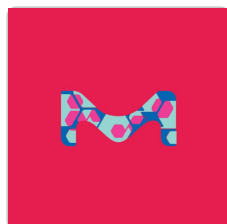
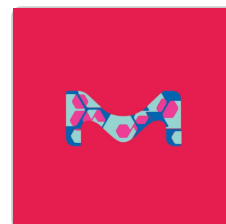
Determination of site-specific glycan heterogeneity on glycoproteins.

Daniel Kolarich et al.*Nature protocols*, 7(7), 1285-1298 (2012-06-09)

The comprehensive analysis of protein glycosylation is a major requirement for understanding glycoprotein function in biological systems, and is a prerequisite for producing recombinant glycoprotein therapeutics. This protocol describes workflows for the characterization of glycopeptides and their site-specific heterogeneity, showing

Comparability analysis of anti-CD20 commercial (rituximab) and RNAi-mediated fucosylated antibodies by two LC-MS approaches.

Chen Li et al.

In developing biosimilar or biobetter products, comparability to the reference product is required to claim similar integrity or intended purpose. In this work, an anti-CD20 monoclonal antibody developed using RNA interference to decrease core fucosylation (RNAi-mediated) was comprehensively characterized by

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Compatible Ends

Enzyme with compatible ends	Recognition sequence	How sequence 5' ends in	How sequence 3' ends in	Enzyme that cuts at this site
BstI	GAATTC	TCTAGC	GAATTC	BstI
XbaI	GAATTC	TCTAGC	GAATTC	XbaI
SmaI	CCCGGG	GGGCCC	CCCGGG	SmaI
EcoRV	GTACCT	GTACCT	GTACCT	EcoRV

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Contactez notre Service technique

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.