[Toutes les photos](#) (1)

Documents

[↓ FDS](#)[🔍 COO/COA](#)[Plus de documents](#) »

NGLYF-RO Roche

N-Glycosidase F

PNGase F of *Flavobacterium meningosepticum*, recombinant from *E. coli*Synonyme(s):
glycosidaseNuméro de classification
(Commission des enzymes): **3.5.1.52** ([BRENDA](#), [IUBMB](#))

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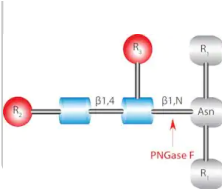
NGLYFL-RO

N-Glycosidase F

recombinant form of the gene from *Flavobacterium meningosepticum*

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Peptide-N-glycosidase F, Peptide-N⁴-(acetyl-β-glucosaminyl)-asparagine amidase

Application

Use N-glycosidase F to cleave all types of asparagine-bound N-glycans, provided that the amino group as well as the carboxyl group are present in a peptide linkage, and that the oligosaccharide has the minimum length of the chitobiose core unit. The reaction products are ammonia, aspartic acid (in the peptide chain), and the complete oligosaccharide.

Note: N-Glycosidase F, recombinant is also available as a lyophilizate without glycerol.

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One unit is the enzyme activity which hydrolyzes 1 nmol dabsyl fibrin glycopeptide or 0.2 nmol dansyl fetuin glycopeptide within 1 minute at 37 °C at pH 7.8.

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Solution in 50 mM sodium phosphate, 12.5 mM EDTA, 50% glycerol (v/v), pH 7.2

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For life science research only. Not for use in diagnostic procedures.

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The sale of the Product does not exhaust or grant any rights in third party patents including patents of companies of the F. Hoffmann - La Roche AG group of companies, in particular, for the use of modified antibodies obtained by using the product.

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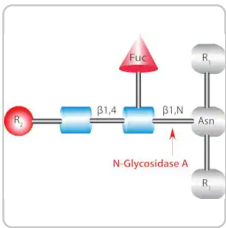
G0535

Glycopeptidase A from almonds

buffered aqueous glycerol solution, ≥0.05 unit/mL

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The table provides a key to monosaccharide symbols used in the product information. It lists various monosaccharides and their corresponding symbols, along with the linkages between them. The symbols include circles, squares, triangles, and diamonds, each representing a different monosaccharide. The linkages are indicated by the orientation of the symbols and the lines connecting them.

Key to Monosaccharide Symbols	
β -D-Glucose (Glc)	β -D-Galactose (Gal)
β -D-Mannose (Man)	α -D-Glucose (Glc) and β -D-Galactose (Gal)
β -D-Galactose (Gal)	β -D-Glucose (Glc) and β -D-Galactose (Gal)
β -D-Mannose (Man)	α -D-Glucose (Glc) and β -D-Galactose (Gal)
β -D-Glucose (Glc)	α -D-Glucose (Glc) and β -D-Galactose (Gal)

Sigma-Aldrich

G1163

O-Glycosidase from *Streptococcus pneumoniae*

recombinant, expressed in *E. coli*, buffered aqueous solution

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Small scale affinity purification and high sensitivity reversed phase nanoLC-MS N-glycan characterization of mAbs and fusion proteins.

Fabian Higel et al.*mAbs*, 6(4), 894-903 (2014-05-23)

N-glycosylation is a complex post-translational modification with potential effects on the efficacy and safety of therapeutic proteins and known influence on the effector function of biopharmaceutical monoclonal antibodies (mAbs). Comprehensive characterization of N-glycosylation is therefore important in biopharmaceutical development. In

New capillary gel electrophoresis method for fast and accurate identification and quantification of multiple viral proteins in influenza vaccines.

Ewoud van Tricht et al.*Talanta*, 144, 1030-1035 (2015-10-11)

Current methods for the identification and/or quantification of viral proteins in influenza virus and virosome samples suffer from long analysis times, limited protein coverage and/or low accuracy and precision. We studied and optimized capillary gel electrophoresis (CGE) in order to

A comparison of anti-HER2 IgA and IgG1 in vivo efficacy is facilitated by high N-glycan sialylation of the IgA.

Gerard Ja Rouwendal et al.*mAbs*, 8(1), 74-86 (2015-10-07)

Monomeric IgA has been proposed as an alternative antibody format for cancer therapy. Here, we present our studies on the production, purification and functional evaluation of anti-HER2 IgA antibodies as anti-cancer agents in comparison to the anti-HER2 IgG1 trastuzumab. MALDI-TOF

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Protocoles

N-Glycosidase F Protocol & Troubleshooting

N-Glycosidase F Protocol & Troubleshooting

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Roche

11363743001**Nutridoma™-CS**

suitable for hybridoma, solution, pkg of 10 mL (50x)

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.