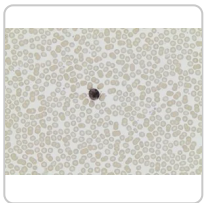
Want a pin for your lab coat?
Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

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91A ▶ Sigma-Aldrich.

α -Naphthyl Acetate (Non-Specific Esterase)

★★★★★ (0)

Kit formulated with all liquid reagents.

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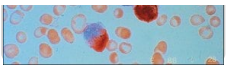
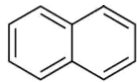
[Cytology Stains](#)

Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

GET YOURS HERE

DESCRIPTION

Application

Esterases are ubiquitous in nature and encompass a series of different enzymes acting upon select substrates. According to the methodology, α -naphthyl acetate is enzymatically hydrolyzed, liberating a free naphthol compound. This then couples with a diazonium compound, forming black colored deposits at sites of non-specific esterase activity. This enzyme is detected primarily in monocytes, macrophages and histocytes, and is normally absent in granulocytes. Lymphocytes may occasionally exhibit enzyme activity. A sodium fluoride inhibition procedure may be run to demonstrate monocyte inhibition. The procedure is run at 37 °C and may be utilized with blood, bone marrow films, tissue-touch preparations, cytocentrifuge preparations and frozen sections. The enzyme does not survive paraffin processing.

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Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

GET YOURS HERE

Reagent product	Description	Download	Price
916	α-Naphthyl acetate solution 10 mL	↓	Tarif
915	Citrate Solution, pH 3.6±0.1 (25 °C), 27 mM 50 mL	↓	Tarif
917	Fast Blue BB Base solution 10 mL	↓	Tarif
914	Sodium nitrite solution 10 mL	↓	Tarif

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916

α-Naphthyl acetate solution

[View Pricing](#)

917

Fast Blue BB Base solution

[View Pricing](#)

914

Sodium nitrite solution

Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

[GET YOURS HERE](#)**915**Citrate Solution, pH 3.6 $\pm$ 0.1 (25 °C), 27 mM[View Pricing](#)**91C**

Naphthol AS-D Chloroacetate (Specific Esterase) Kit

[View Pricing](#)**Produit Comparable****90A1** α -Naphthyl Acetate (Non-Specific Esterase), Select reagents packaged in gelatin capsules.[View Pricing](#)**INFORMATIONS SUR LA SÉCURITÉ****Pictograms****GHS02,GHS05,GHS06,GHS08****Mention d'avertissement**

Danger

Mentions de danger**H226 - H290 - H301 - H312 +
H332 - H319 - H370 - H412****Conseils de prudence****P210 - P273 - P280 - P301 +
P310 - P303 + P361 + P353 -
P304 + P340 + P312**

Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

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e.g. 023J5431

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Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

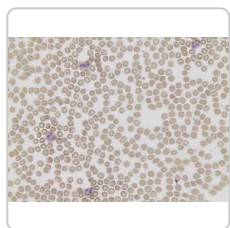
GET YOURS HERE

Plus De Documents

NAPHTHOL AS-D CHLOROACETATE ESTERASE AND α -NAPHTHYL ACETATE ESTERASE FDS

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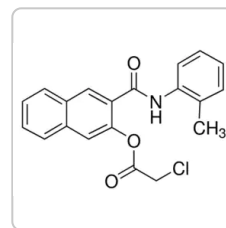
Sigma-Aldrich

387A

Acid Phosphatase, Leukocyte (TRAP) Kit
Kit formulated with all liquid reagents

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SA Kuznetsov et al.

The Journal of cell biology, 153(5), 1133-1140 (2001-05-31)

We report the isolation of adherent, clonogenic, fibroblast-like cells with osteogenic and adipogenic potential from the blood of four mammalian species. These cells phenotypically resemble but are distinguishable from skeletal stem cells found in bone marrow (stromal stem cells, "mesenchymal

Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

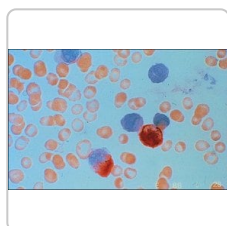
GET YOURS HERE

The Journal of investigative dermatology, 135(1), 279-288 (2014-07-23)

During wound healing, excessive inflammation, angiogenesis, and differentiated human dermal fibroblast (HDF) function contribute to scarring, whereas hyperpigmentation negatively affects scar quality. Over 100 million patients heal with a scar every year. To investigate the role of the beta 2

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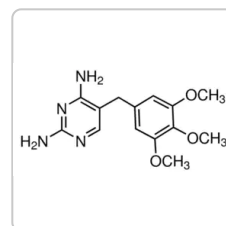
Contactez notre Service technique

Want a pin for your lab coat?

Subscribe to our scientific newsletters to get a shiny 'Magical Flask' pin.

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.