Brighten up your lab coat!

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Description générale

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

of novel microbial species. Applications of metagenomic studies include public health data analysis, discovery of novel proteins, enzymes and natural products, environmental studies, and agricultural investigations.

Application

MetaPolyzyme is a mixture of 6 enzymes that is intended for isolation of total DNA for metagenomics studies. Adapted from an initial formulation devised by S. Tighe,² MetaPolyzyme was evaluated and developed in consultation and collaboration with the Association of Biomolecular Resource Facilities (ABRF) Metagenomics and Microbiome Research Group (MMRG). Test data on the use of MetaPolyzyme is accessible at the [2017 ABRF MGRG Poster](#).

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Microbes can be difficult to disrupt because the cell walls may form capsules or resistant spores. DNA can be extracted by using lysing enzymes (such as lyticase or chitinase) to induce partial spheroplast formation. Spheroplasts are subsequently lysed to release DNA. MetaPolyzyme is intended for use in metagenomic studies, such as study of microbes and microbiomes of extreme and unique environments. MetaPolyzyme is intended for digestion of microbes for evaluation by whole genome shotgun sequencing for metagenomics and metatranscriptomic approaches.

Composants

The enzymes in MetaPolyzyme are:

- Achromopeptidase
- Chitinase
- Lyticase
- Lysostaphin
- Lysozyme

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

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e.g. 023J5431

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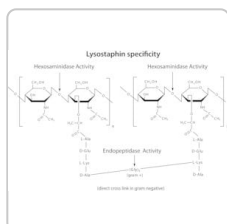
Rechercher

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Product Information Sheet - MAC4L FDS

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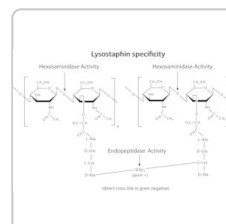
L4402

Lysostaphin from *Staphylococcus staphylolyticus*

BioUltra, ≥97% (SDS-PAGE), Protein 40-60 % by biuret, ≥2,000 units/mg protein

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SRE0053

Lysostaphin from *Staphylococcus staphylolyticus*

Suitable for manufacturing of diagnostic kits and reagents., lyophilized powder

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Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

Microbiomic Analysis on Low Abundant Respiratory Biomass Samples; Improved Recovery of Microbial DNA From Bronchoalveolar Lavage Fluid.

Saladie M et al.

Frontiers in Microbiology, 11, 572504-572504 (2020)

Erratum to: The Metagenomics and Metadesign of the Subways and Urban Biomes (MetaSUB) International Consortium Inaugural Meeting Report.

Microbiome, 4(1), 45-45 (2016-08-20)

Metagenome sequencing-based strain-level and functional characterization of supragingival microbiome associated with dental caries in children.

Nezar Noor Al-Hebshi et al.

Journal of oral microbiology, 11(1), 1557986-1557986 (2019-01-24)

Studies of the microbiome associated with dental caries have largely relied on 16S rRNA sequence analysis, which is associated with PCR biases, low taxonomic resolution, and inability to accurately study functions. Here, we employed whole metagenome shotgun sequencing, coupled with

Cartography of Opportunistic Pathogens and Antibiotic Resistance Genes in a Tertiary Hospital Environment.

Chng KR et al.

Nature Medicine, 26(6), 941-951 (2020)

Genomic Methods and Microbiological Technologies for Profiling Novel and Extreme Environments for the Extreme Microbiome Project (XMP).

Scott Tighe et al.

Journal of biomolecular techniques : JBT, 28(1), 31-39 (2017-03-25)

The Extreme Microbiome Project (XMP) is a project launched by the Association of Biomolecular Resource Facilities Metagenomics Research Group (ABRF MGRG) that focuses on whole genome shotgun sequencing of extreme and unique environments using a wide variety of biomolecular techniques.

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

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Articles

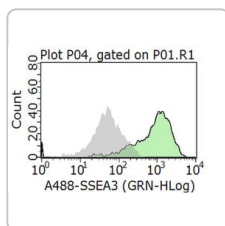
Lysing Enzymes

Enzymes provide a non-mechanical method for cell lysis and protoplast preparation. It may seem like a simple process to throw in your enzyme, stick your tube in the water-bath and walk away, but what is actually going on in that process?

Human Microbiome Research Workflow: An Overview of the Role in Human Health and Disease

An overview of human microbiome research, workflow challenges, sequencing, library production, data analysis, and available microbiome reagents to support your research.

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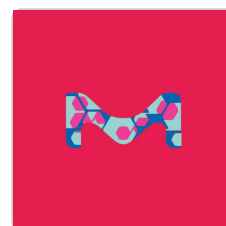
Sigma-Aldrich

MAB4303-I

Anti-Stage-Specific Embryonic Antigen-3
Antibody, clone MC-631
clone MC-631, from rat

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Sigma-Aldrich

M9956

M9 Minimal Salts, 5X
Liquid microbial growth medium

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.