Brighten up your lab coat!

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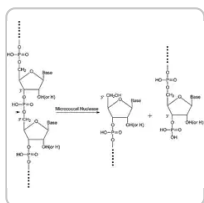
🔍 COO/COA

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N3755 ▶ **Sigma-Aldrich**

Nuclease micrococcal from *Staphylococcus aureus*

★★★★★ (0)

100-300 units/mg protein

Synonyme(s):

Micrococcal nuclease, Endonuclease micrococcal, MNase, Micrococcal endonuclease

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[Ajouter au panier](#)

Brighten up your lab coat!

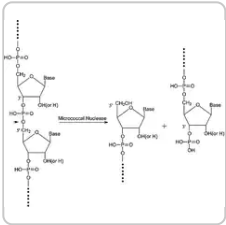
Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

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
Sigma-Aldrich
N5386
Nuclease micrococcal from *Staphylococcus aureus*
100-300 units/mg protein

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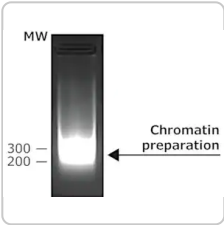Sigma-Aldrich
CS0004
Nuclease micrococcal from *Staphylococcus aureus*
With reaction buffer and BSA solution

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DESCRIPTION

Application

Nuclease from *Staphylococcus aureus* has been used in a study to assess coagulase and heat-resistant strains found in animals. It has also been used in a study to investigate the expression characteristic of two genes in *S. aureus* that encode two thermostable nucleases.

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Cleavage of DNA by Nuclease micrococcal is greater at the 5' sides of A or T nucleotides than at G or C.

Hydrolyzes 5'-phosphodiester bonds of DNA.

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Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

[GET YOURS HERE](#)**Autres remarques**

Note: One μ molar unit = approx. 85 A_{260} units, where an A_{260} unit is equivalent to a ΔA_{260} of 1.0 in 30 min at pH 8.8 at 37 °C in a 3.55 mL reaction volume. Light path = 1 cm.

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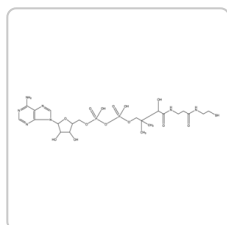
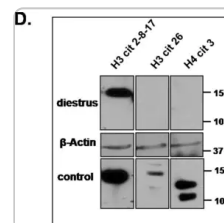
Sigma-Aldrich

07-596

Anti-Histone H4 (citruiline 3) Antibody

Upstate®, from rabbit

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

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Comparative expression analysis of two thermostable nuclease genes in *Staphylococcus aureus*.

Yu Hu et al.

Foodborne pathogens and disease, 9(3), 265-271 (2012-03-03)

Thermonuclease is known as a specific virulence factor in *Staphylococcus aureus*. It is widely used as a genetic marker for detection of *S. aureus* in various types of food. Previous studies have revealed the existence of two functional thermostable nucleases

High sequence specificity of micrococcal nuclease.

C Dingwall et al.

Nucleic acids research, 9(12), 2659-2673 (1981-06-25)

The substrate specificity of micrococcal nuclease (EC 3.1.4.7.) has been studied. The enzyme recognises features of nucleotide composition, nucleotide sequence and tertiary structure of DNA. Kinetic analysis indicates that the rate of cleavage is 30 times greater at the 5'

Coagulase and Heat-resistant Nuclease Producing *Staphylococcus epidermidis* Strains from Animals

DEVRIESE, L. and P. OEDING

Journal of Applied Microbiology, 39, 197-207 (1975)

Noninvasive imaging of *Staphylococcus aureus* infections with a nuclease-activated probe.

Frank J Hernandez et al.

Nature medicine, 20(3), 301-306 (2014-02-04)

Technologies that enable the rapid detection and localization of bacterial infections in living animals could address an unmet need for infectious disease diagnostics. We describe a molecular imaging approach for the specific, noninvasive detection of *S. aureus* based on the

Intracellular nucleosomes constrain a DNA linking number difference of -1.26 that reconciles the Lk paradox.

Joana Segura et al.

Nature communications, 9(1), 3989-3989 (2018-09-30)

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

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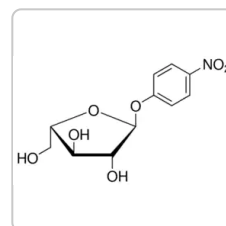
Sigma-Aldrich

N3641

4-Nitrophenyl α-L-arabinofuranoside

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