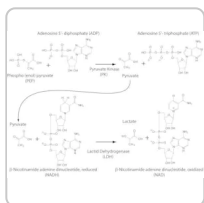
Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

[GET YOURS HERE](#)

Toutes les
photos (1)

Documents

FDS

COO/COA

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P0294 ▶ **Sigma-Aldrich.**

Pyruvate Kinase/Lactic Dehydrogenase enzymes from rabbit muscle

★★★★★ (0)

For the Determination of ADP, buffered aqueous glycerol solution

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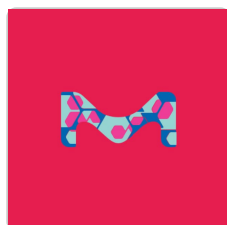
Brighten up your lab coat!

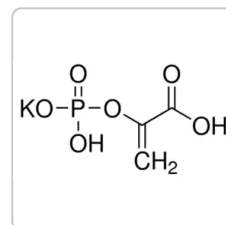
Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

[GET YOURS HERE](#)

Sigma-Aldrich

P1506**Pyruvate Kinase from rabbit muscle**

Type II, ammonium sulfate suspension, 350-600 units/mg protein

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Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

Description générale

Pyruvate Kinase from rabbit muscle is a metalloenzyme which catalyzes the conversion of phosphoenol pyruvate to pyruvate in the glycolysis pathway. It corresponds to a molecular weight of 59 kDa. It exists as a tetramer and undergoes conformational changes in the active site to accommodate substrate. Lactic dehydrogenase (LDH) catalyzes the lactate to pyruvate conversion in anaerobic glycolysis. It exists as tetramer and comprises of two subunits (H and M). The LDH of eukaryotes undergo active-site loop gating for their catalytic functionality.

Application

Pyruvate Kinase/Lactic Dehydrogenase enzymes from rabbit muscle has been used:

- for ATP generation in the active microtubule preparation
- in the enzyme linked ATPase assay of skeletal muscle heavy meromyosin (HMM)
- as a standard control for quantifying mesenchymal stem cells (MSCs) lactate dehydrogenase

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ADP Quantification Assay protocol for the use of PK/LDH in the determination of ADP. Solutions containing unknown concentrations of ADP can be substituted for reagent D in this protocol. Further dilutions of the ADP solution may be required

Lactate dehydrogenase from rabbit muscle can be inhibited by ascorbate. Aldolase and actin were shown to block this inhibitory effect.

Pyruvate kinase also catalyzes the phosphorylation of thiamine diphosphate (TDP) to thiamine triphosphate (TTP) which may find application in antiviral and tumor therapy.

Pyruvate kinase requires bivalent and monovalent cations such as Mg^{2+} and K^{+} respectively for activation to occur.

Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

[GET YOURS HERE](#)**Forme physique**

Solution in 50% glycerol containing 10 mM HEPES, pH 7.0, 100 mM KCl and 0.1 mM EDTA

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Eyeshields, Gloves, multi-purpose combination respirator cartridge (US)

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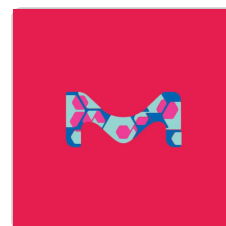
Sigma-Aldrich

P1903

Pyruvate Kinase from *Bacillus stearothermophilus*

Type VIII, lyophilized powder, 100-300 units/mg protein

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Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

Targeting Bcr-Abl by combining allosteric with ATP-binding-site inhibitors.

Jianming Zhang et al.

Nature, 463(7280), 501-506 (2010-01-15)

In an effort to find new pharmacological modalities to overcome resistance to ATP-binding-site inhibitors of Bcr-Abl, we recently reported the discovery of GNF-2, a selective allosteric Bcr-Abl inhibitor. Here, using solution NMR, X-ray crystallography, mutagenesis and hydrogen exchange mass spectrometry

The T790M mutation in EGFR kinase causes drug resistance by increasing the affinity for ATP.

Cai-Hong Yun et al.

Proceedings of the National Academy of Sciences of the United States of America, 105(6), 2070-2075 (2008-01-30)

Lung cancers caused by activating mutations in the epidermal growth factor receptor (EGFR) are initially responsive to small molecule tyrosine kinase inhibitors (TKIs), but the efficacy of these agents is often limited because of the emergence of drug resistance conferred

Ligand-Induced Domain Movement in Pyruvate Kinase: Structure of the Enzyme from Rabbit Muscle with Mg²⁺, K⁺, and I-Phospholactate at 2.7 Å Resolution

Larsen TM, et al.

Archives of Biochemistry and Biophysics, 345(2), 199-206 (1997)

Thermal activation of `allosteric-like?large-scale motions in a eukaryotic Lactate Dehydrogenase

Katava M, et al.

Scientific Reports, 7, 41092-41092 (2017)

Chemical and enzymatic characterization of recombinant rabbit muscle pyruvate kinase

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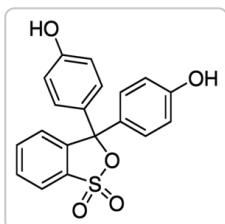
Brighten up your lab coat!

Subscribe to our scientific newsletters and make your lab mates jealous with a shiny 'Magical Flask' pin.

GET YOURS HERE

Sigma-Aldrich

P0290

Phenol red solution

0.5%, liquid, sterile-filtered, BioReagent, suitable for cell culture

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.