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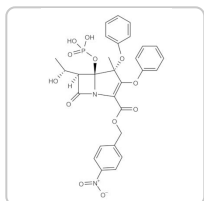
COO/COA

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P6556 ► Sigma-Aldrich.

Proteinase K from *Tritirachium album*

★★★★★ (0)

lyophilized powder, ≥30 units/mg protein

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Sigma-Aldrich

P2308**Proteinase K from *Tritirachium album***

lyophilized powder, BioUltra, ≥30 units/mg protein, for molecular biology

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Millipore

70663**Proteinase K, Lyophilized**

Highly active serine protease that exhibits broad cleavage specificity on native and denatured proteins...

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DESCRIPTION

Application

Product P6556 is provided as a lyophilized powder. Product P6556 has been used to break down human lens protein. Protease footprinting by Proteinase K digestion can reveal protein-protein surface interactions. The enzyme from Sigma has been used in the pre-hybridization step of chicken embryos. It has also been used for the enrichment of PrP^{Sc}, a prion protein that is present in sheep, hamster and mouse scrapie samples.

Proteinase K is useful for the proteolytic inactivation of nucleases during the isolation of DNA and RNA. It is used for the removal of endotoxins bound to cationic proteins such as lysozyme and ribonuclease A. It is useful for the isolation of hepatic, yeast, and mung bean mitochondria

and is used to determine enzyme localization on membranes

It is used for the treatment of paraffin embedded tissue sections to expose antigen binding sites for antibody labeling and for digestion of proteins from brain tissue samples for prions in Transmissible Spongiform Encephalopathies (TSE) research. Product P6556 is provided as a lyophilized powder. Product P6556 has been used to break down human lens protein.

Useful for the proteolytic inactivation of nucleases during the isolation of DNA and RNA.

Removes endotoxins that bind to cationic proteins such as lysozyme and ribonuclease A.

Reported useful for the isolation of hepatic, yeast, and mung bean mitochondria

Determination of enzyme localization on membranes

Treatment of paraffin embedded tissue sections to expose antigen binding sites for antibody labeling.

Digestion of proteins from brain tissue samples for prions in Transmissible Spongiform Encephalopathies (TSE) research.

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Proteinase K has a broad specificity and degrades many proteins even in the native state. It mainly cleaves the peptide bond adjacent to the carboxyl group of aliphatic and aromatic amino acids with blocked α -amino groups. The optimum pH is between 7.5-9.0 and the isoelectric point is 8.9 Ca^{2+} (1-5 mM) is required for activation. Proteinase K is inhibited by diisopropyl fluorophosphate (DFIP), and phenylmethanesulfonyl fluoride (PMSF).

Proteinase K is a stable and highly reactive serine protease. Evidence from crystal and molecular structure studies indicates the enzyme belongs to the subtilisin family with an active-site catalytic triad (Asp^{39} - His^{69} - Ser^{224}). It is stable in a broad range of environments: pH, buffer salts, detergents (SDS), and temperature. In the presence of 0.1-0.5% SDS, proteinase K retains activity and will digest a variety of proteins and nucleases in DNA preparations without compromising the integrity of the isolated DNA.

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One unit will hydrolyze urea-denatured hemoglobin to produce color equivalent to 1.0 μ mole of tyrosine per min at pH 7.5 at 37 °C (color by Folin-Ciocalteu reagent).

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dust mask type N95 (US),
Eyeshields, Faceshields,
Gloves

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
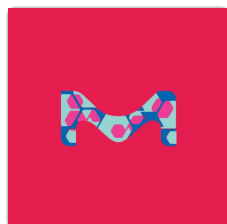
1.07393

Proteinase K

(from *Tritirachium album*) solution in Tris/HCl pH 7.5;
0.01 mol/l; 600 mAnson-U/ml; for molecular biology EC...

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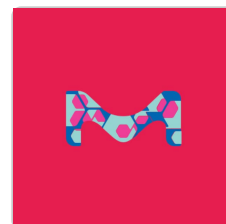Structure elucidation of a novel yellow chromophore from human lens protein.

Rongzhu Cheng et al.

The Journal of biological chemistry, 279(44), 45441-45449 (2004-08-19)

We report here the isolation of a novel acid-labile yellow chromophore from the enzymatic digest of human lens proteins and the identification of its chemical structure by liquid chromatography-mass spectrometry, liquid chromatography-tandem mass spectrometry, and (1)H, (13)C, and two-dimensional NMR.

Properties of a subtilisin-like proteinase from a psychrotrophic *Vibrio* species comparison with proteinase K and aqualysin I.

M M Kristjánsson et al.

European journal of biochemistry, 260(3), 752-760 (1999-04-02)

An extracellular serine proteinase purified from cultures of a psychrotrophic *Vibrio* species (strain PA-44) belongs to the proteinase K family of the superfamily of subtilisin-like proteinases. The enzyme is secreted as a 47-kDa protein, but under mild heat treatment (30

Stimulation of proteinase K action by denaturing agents: application to the isolation of nucleic acids and the degradation of 'masked' proteins.

H Hilz et al.

European journal of biochemistry, 56(1), 103-108 (1975-08-01)

Hydrolysis of serum albumin by proteinase K was strongly (greater than 7-fold) stimulated by urea and dodecylsulfate in a dose-dependent manner. With an oligopeptide as substrate, however, proteinase K was inactivated by dodecylsulfate. This indicates that the apparent activation of

FISHing for chick genes: Triple-label whole-mount fluorescence in situ hybridization detects simultaneous and overlapping gene expression in avian embryos.

Nathaniel Denkers et al.

Developmental dynamics : an official publication of the American Association of Anatomists, 229(3), 651-657 (2004-03-03)

Multi-color whole-mount in situ hybridization is a powerful technique for comparing the spatial expression patterns of two or more genes in developing embryos. We have developed an amplified triple-label whole-mount fluorescence in situ hybridization (FISH) protocol that permits detection of

Gonadotropin characterization, localization and expression in the European hake (*Merluccius merluccius*).

Michela Candelma et al.

Reproduction (Cambridge, England), 153(2), 123-132 (2016-11-03)

In vertebrates, the regulation of gametogenesis is under the control of gonadotropins (Gth), follicle-stimulating hormone (Fsh) and luteinizing hormone (Lh). In fish, the physiological role of Gths is not fully understood, especially in species with asynchronous ovarian development. To elucidate

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Articles

Enzymatic Assay of Proteinase K with Hemoglobin Substrate

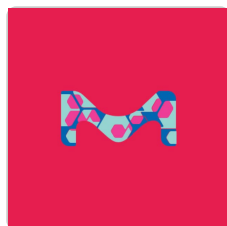
Proteinase K (EC 3.4.21.64) activity can be measured spectrophotometrically using hemoglobin as the substrate. Proteinase K hydrolyzes hemoglobin denatured with urea, and liberates Folin-positive amino acids and peptides. One unit will hydrolyze hemoglobin to produce color...

Protocoles

Enzymatic Assay of Proteinase K (EC 3.4.21.64)

Proteinase K (EC 3.4.21.64) activity can be measured spectrophotometrically using hemoglobin as the substrate. Proteinase K hydrolyzes hemoglobin denatured with urea, and liberates Folin-positive amino acids and peptides. One unit will hydrolyze hemoglobin to produce color...

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
Sigma-Aldrich

P6534

Phospholipase A2 from porcine pancreas
ammonium sulfate suspension, ≥600 units/mg protein

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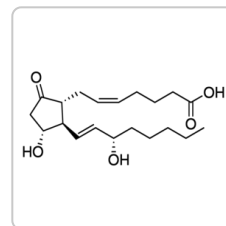Sigma-Aldrich

P6532

Prostaglandin E2
γ-irradiated, powder, BioXtra, suitable for cell culture

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Contactez notre Service technique

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© 2022 Merck KGaA, Darmstadt, Allemagne et/ou ses sociétés affiliées. Tous droits réservés.